



## Quick - Start

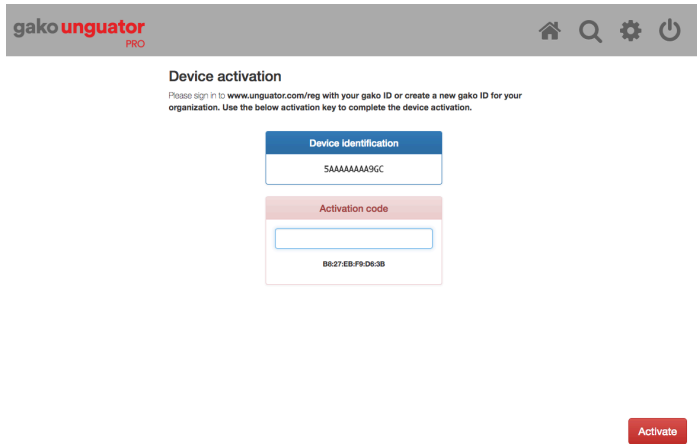

### How to register your gako unguator PRO device

- Go to [unguator.com/reg](http://unguator.com/reg)
- Enter the 12 digit serial number
- Enter the activation code on your device

--> review chapter 3

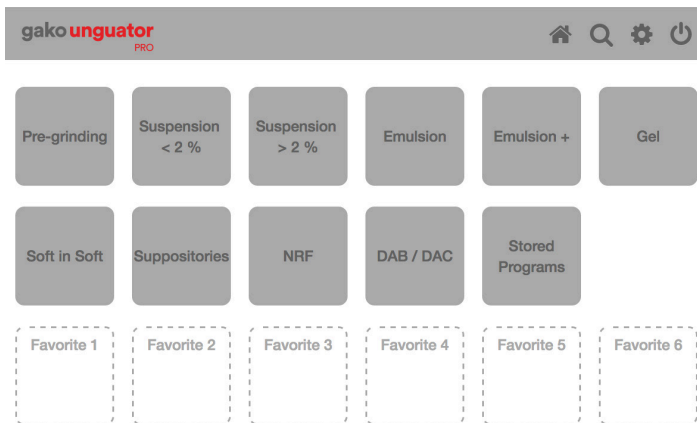


### Display handling

- Touch screen reacting to pressing directly the respective symbols with your finger
- main display displaying the mixing programs
- after selecting the program, scan of the jar size via the integrated camera or manual input by selecting „jar size“

--> review chapter 3

### Lift arm with quick-lock coupling

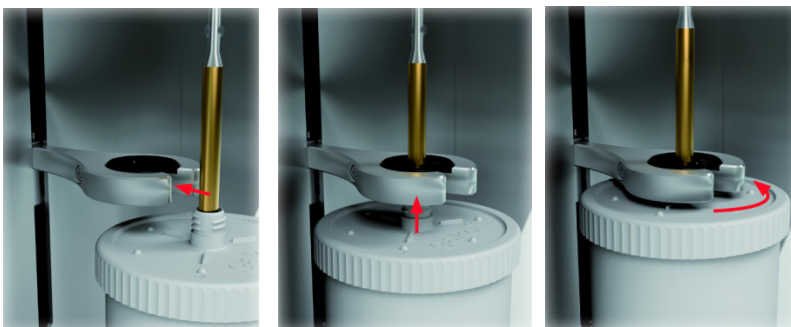

- Hold the gako logo on top of the jar to the front; insert the mixing shaft directly from the front into the lift arm
- Insert the mixing unit from below
- Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lift arm and push the start button

--> review chapter 4