



## Quick - Start

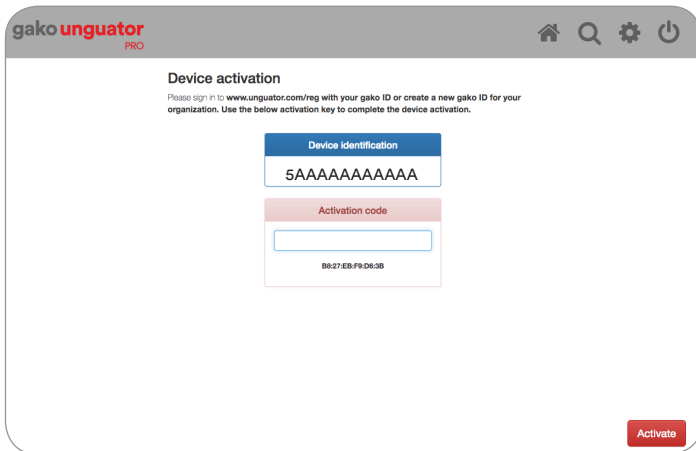

### How to register your gako unguator PRO device

Go to [unguator.com/reg](http://unguator.com/reg)  
Enter the 12 digit serial number  
Enter the activation code on your device

--> review chapter 3

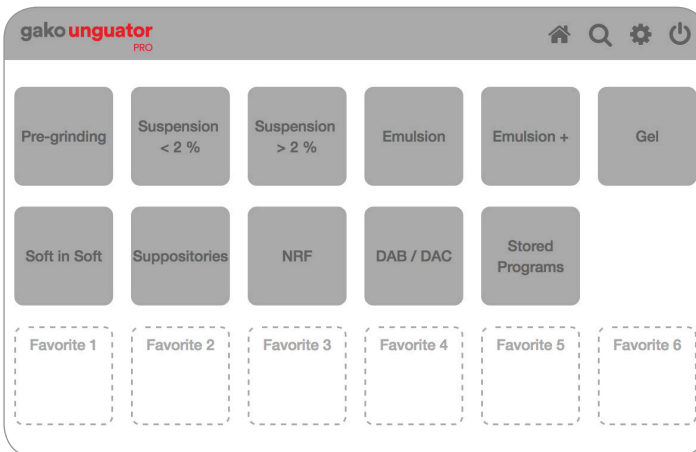


### Display handling

Touch screen reacting to pressing directly the respective symbols with your finger  
Main display displaying the mixing programs  
After selecting the program, scan of the jar size via the integrated camera or manual input by selecting „jar size“

--> review chapter 3

### Lifting arm with quick-lock coupling

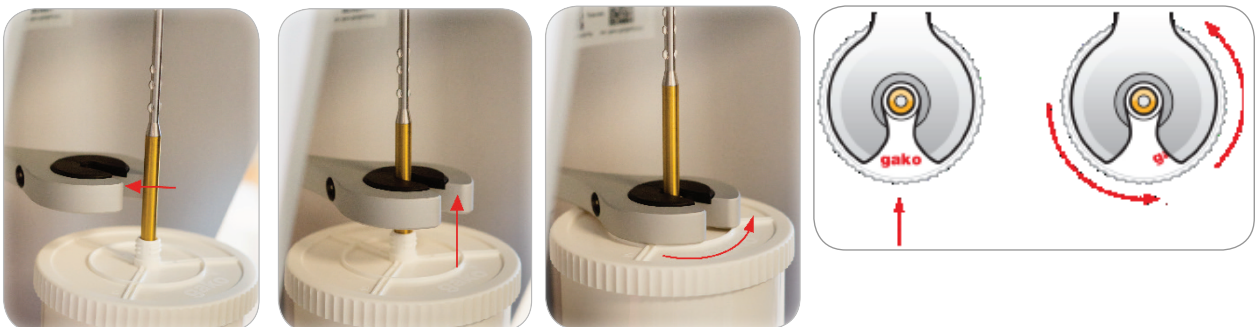

- Hold the gako logo on top of the jar to the front; insert the mixing shaft directly from the front into the lifting arm
- Insert the mixing unit from below
- Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lifting arm and push the start button

--> review chapter 4