
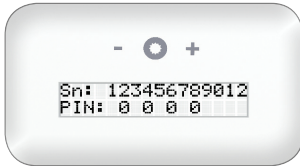



Quick - Start

How to register your gako unguator BASIC device

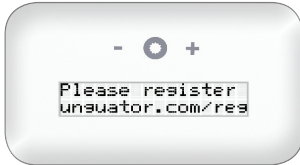


- Go to **unguator.com/reg**
- Enter the 12 digit serial number
- Enter the four digit unlock PIN on your device
--> review chapter 3

Display handling

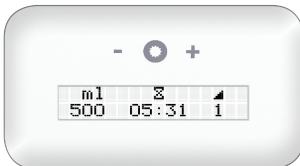

- Touch panel with minus button (left), gako button (middle), plus button (right)
- start display shows jar size, mixing time and mixing speed
- After pushing the gako button for two seconds “Start” appears on the display and the jar can be mounted
--> review chapter 3

Lifting arm with integrated quick-lock coupling

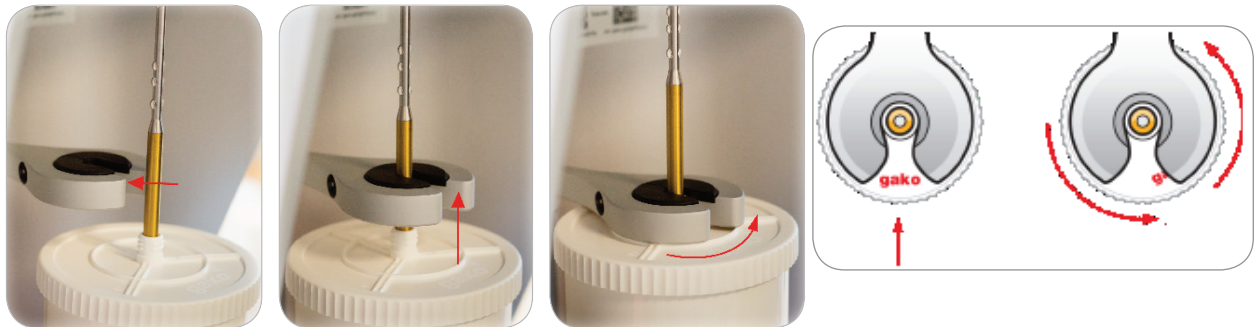

- Hold the gako logo on top of the jar to the front; insert the mixing shaft directly from the front into the lifting arm
- Insert the mixing unit from below
- Apply slight pressure from the top onto the lifting arm. Fix the jar with a slight turn counterclockwise (about a quarter turn) into the lifting arm
- Move the mixing unit in the lifting arm by hand all the way up and push the gako button
--> review chapter 3

Video link, how to mount the jar
goo.gl/gtDmLt

